



**HISTOPen**

Catalog #: 30020

[www.histopat.hu](http://www.histopat.hu)

**DATA SHEET**

<b>Description:</b>	HISTOPen provides a thin, film-like water repellent barrier, when a circle is drawn around the specimen on a glass slide. This barrier creates a proper surface tension to hold the applied reagents within the target area on the slide. Two specimen separated by two circles can be applied to the same slide.
<b>Intended use:</b>	For histology, cytology, immunoenzymatic-, immunofluorescent-, ISH-methods, microbiological- and virological tests.
<b>Properties:</b>	<ol style="list-style-type: none"><li>1. Heat resistant up to 120 °C</li><li>2. Ready to use on slides</li><li>3. Buffer (water)-resistant</li><li>4. Alcohol-resistant</li><li>5. Soluble in xylene!</li></ol>
<b>Application:</b>	Before the first use, push the pen gently until the liquid saturates the tip properly. Before the application of the antibody draw a circle around the sample on the glass slide.
<b>Storage:</b>	Store with tightly closed cap in horizontal position.

**For Research Use Only!**

